

## **KL-03(II)A PH Meter, Lab China Economical Cheaper PH Teter**

### **Key Features:**

Pocket sized and feather-weight for portable use in a large number of applications. It is utilize user-replaceable electrodes.

### **Specification:**

Measuring range: 0.00-14.00pH

Accuracy:  $\pm 0.1$ PH

Resolution: 0.01PH

Operating temperature : 0C- 50C

Automatic temperature compensation: 0C- 50C

Plug: BNC+1m cable

Dimensions: 120mm x 26mm x 26mm

Weight: 70g

lab china economical cheaper ph meter

### **Operation**

- 1.Remove the protective cap of electrode.
- 2.First rinse the electrode with distilled water, and suck it with filter-paper.
- 3.Turn the meter on by pressing the ON key.
- 4.Immerse the ph meter electrode in the solution to be tested.
- 5.Stir gently and wait for the reading to stabilize.
- 6.After use, rinse the electrode with water to minimize contamination. Turn the meter on by pressing the OFF key.
- 7.Always replace the protective cap after use.

lab china economical cheaper ph meter