

PH-01B Promotional Digital Pen Type PH Meter Manufacturer

Specification:

1. Measuring ranges:0.0-14.0pH
2. Accuracy:±0.01PH
3. Resolution: 0.01PH
4. Operating temperature : 0~80°C
5. Calibration: two points automatic calibration
6. Power supply:2*1.5v(keep using more than 500 hours)
7. Dimension: 153*29*16mm
8. Weight: 52g

Operation:

- 1.Remove the protective cap.
- 2.First rinse the electrode with distilled water ,and dry it with filter pater.
- 3.Turn the meter on by pressing the ON/OFF key.
- 4.Immerse the ph meter electricle in the solution to be tested .
5. Stir gently and wait for the reading to stablilize.
- 6.After finished ,clear the electrode with distilled water turn the meter off by pressing the "ON/OFF" key.
- 7.Always replace the protective cap after use .